🛟 eurofins

GeneScan

🛟 eurofins

GeneScan

DN*Animal Ident* Sheep

IPC (LR/HR)

Cat. No. 5422211710

Test kit for the qualitative detection of sheep DNA with IPC, 96 real-time PCR reactions

DNAnimal Ident Sheep IPC (LR/HR)_ID2554

V3 19.01.2018

🛟 eurofins

GeneScan

Table of Contents:

4
7
8
9
IN g
10
12
12
13
15
17

13. TECHNICAL SERVICE

If you have any questions or experience any difficulties regarding this kit or Eurofins GeneScan Technologies products in general, please do not hesitate to contact us. Eurofins GeneScan customers are a major source of information regarding advanced or specialised uses of our products. This information is helpful to other scientists as well as to the researchers at GeneScan Technologies. We therefore encourage you to contact us if you have any suggestions about product performance or new applications and techniques. For technical assistance and more information please call the Eurofins GeneScan Technologies Technical Service Department or your local distributor.

Eurofins GeneScan Technologies GmbH

Engesser Str. 4 79108 Freiburg, Germany Phone: + 49-(0)761-5038-100 Fax: + 49-(0)761-5038-111 kits@eurofins.com www.eurofins.com/kits

© 2018 Eurofins GeneScan Technologies GmbH, all rights reserved.

DNAnimal Ident

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

8. RESULTS	21
8.1 Evaluation	21
8.2 Interpretation of Results	23
8.3 Ambiguous results	27
8.4 Limit of detection	28
9. PRODUCT USE LIMITATIONS	28
10. PRODUCT WARRANTIES AND SATISFACTION GUARANTEE	29
11. IMPORTANT NOTES	30
12. TROUBLESHOOTING	31
13. TECHNICAL SERVICE	32

Test kits, their components and instructions for use are subject to alterations. They are intended for research purposes only.

DNAnimal Ident

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

DNAnimal Ident Sheep IPC Kit

Kit for qualitative real-time PCR detection of sheep DNA

1. INTRODUCTION

National regulations in most countries of the world require declaration of components of food and feed and a corresponding labelling.

Declaration of components of animal origin are crucial e.g. for vegetarian food, food for religious communities, general authenticity and quality control as well as for export and trade.

The Eurofins GeneScan DNAnimal kits comprise several major advantages:

- High sensitivity (LOD 0.01%)
- High specificity (primers and probe)
- · Robust test methods
- Fast results

4/32

🔅 eurofins

GeneScan

The DNAnimal Ident Sheep IPC kit was designed for a general detection of sheep (Ovis aries) DNA in DNA extracted i.e. from food and feed matrices.

The complete test comprises the following steps:

- 1. DNA extraction (not included in this kit)
- 2. Real-time PCR detection of the target

Animal species can be analysed in most matrices, even in gelatine. The processing steps commonly used in food and feed production usually reduce the fragment length of the sample DNA, but DNA is not completely degraded and the base sequence of the fragments is not changed. However, extremely rigid (acid/heat) processing can degrade DNA in a way that amplification is impossible.

Specific DNA sequences are amplified by RT-PCR (Real-time Polymerase Chain Reaction) and detected with high specificity due to the Taqman[™] probes.


The probe of the detection system in this kit uses FAM® as reporter dye and the IPC (internal positive control) uses $R6G^{TM}$. The probes use non-fluorescent quenchers.

The kit is validated for use on Agilent Mx3005P/3000P, ABI7500/Fast, Roche LC480, Bio-Rad CFX96 Touch, AriaMx and ABI 7900HT.

DNAnimal Ident


Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

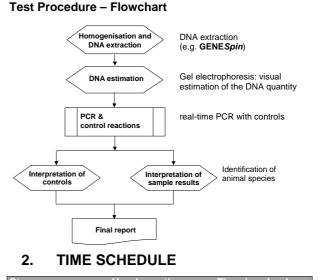

StepHands-on timeTime involvedDNA-extraction0.5 h1.5 hPCR-reaction0.5 h3.5 hTotal time1.0 h5.0 h

The hands-on time has been calculated for a single sample. When several samples are handled simultaneously, the time increases.

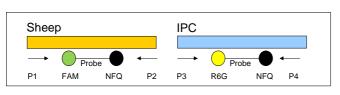

FAMTM and R6G are the fluorescent reporter dyes attached to the 5' ends of the probes for the sheep target and for the IPC (internal positive control). Non-fluorescent quenchers are used for quenching.

Important note:

Please note, cross-reaction with chamois and oryx can occur.

5/32

DNAnimal Ident

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

3. COMPONENTS OF THE KIT

The kit contains all components to run control and specific reactions for a total of 100 PCR reactions.

2x MasterMix QL RT IPC (LR/HR+) GS-P-08.040, Sheep

1 mL composed of

- 650 μL BasicMix QL RT (NR) GS-P-26.012 EFGi TP 2x (NR) 1.5 U
- 390 μL OligoMix QL RT IPC (LR/HR+)

GS-P-08.040 • Sheep

Mix prior to use!

- 2 x Positive control DNA: Genomic DNA from sheep (50 µL, 10 copies/µL)
- 1x DNA stabilization buffer II, 150 µL (for NTCs)
- 2x **ROX™ dye**, 23 μL (to be used only for High Rox cyclers, see instructions below)

Important Note: Store all following components at -20°C. Never store solutions and materials for DNA extraction together with samples or amplicons. Never use them in areas where PCR products (amplicons) may be present.

8 / 32

🛟 eurofins

GeneScan

6. SAMPLE PREPARATION

The extraction of DNA from the sample is a crucial step, and in order to guarantee optimal quantity and purity, an appropriate method has to be chosen for each sample matrix. For most products, the DNA can be extracted with a kit from our DNA*Extractor* kit line or with our GENE*Spin* kit.

Absolute DNA amount and purity (meaning: free of PCR inhibiting substances) affect overall sensitivity of the analysis. Low DNA amount subjected to analysis results in poor LOD with regard to the specific test sample. Insufficiently pure DNA displaying inhibitory effects may even make testing impossible.

It is recommended to use the sample DNA undiluted. In case of inhibitors present in the DNA a repeat of the analysis with diluted sample DNA is feasible, if DNA concentration allows.

4. PREPARATION OF THE KIT'S REAGENTS

Store the kit at -20°C until opened for the first time. Thaw reagents just before use, mix them by vortexing and centrifuge briefly before use. Just thaw as many BasicMix and OligoMix vials as necessary. Frequent freezing and thawing might cause inactivation of the reagents. If smaller volumes are needed, aliquot reagents at first use. Refer to the reagent label for specific instructions regarding the correct storage.

5. MATERIAL AND EQUIPMENT NOT INCLUDED IN THE KIT

- water, DNase-free
- Vortex mixer
- Micropipettes (variable 1-1000 μ L) and filter tips
- PCR optical tubes or plates 0.2 mL and optical caps or seals
- RT-PCR Thermocycler

9 / 32

DNAnimal Ident Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

DNA amount can be determined prior to analysis by one of the following methods:

Real-time PCR monitor run (preferred method)
Fluorimetric or spectrophotometric or other physical measurements, e.g. gel electrophoresis may be applied.

Especially the spectrophotometric methods can suffer significant errors due to interfering DNA impurities or partial DNA denaturation/degradation. For samples known with regard to extraction yield and purity, DNA measurement may be omitted.

Each sample should be treated in duplicate. It is recommended to perform an extraction control for each set of samples extracted simultaneously, i.e. a complete DNA extraction without sample material, which should subsequently undergo PCR analysis.

Cat. No. 5422211710

7. PCR

7.1 Special precautions during PCR

PCR is an exponential reaction. Detection of single target DNA copies is possible. The extreme sensitivity requires special precautions for handling and equipment. A successful amplification leads to several billion amplicons in the reaction tube. Each of them might lead to a false positive result when contaminating sample material or PCR, e.g. by spreading in aerosols. Most important rules to avoid false-positive results are:

a) Separate the different procedures spatially. Use

- a) Separate the different procedures spatially. Use separate rooms for sample preparation, amplification and analysis of amplicons and dedicate separate equipment and materials for each procedure. Mixing of PCR components should be done in a separate room, best on a clean bench. Keep all working steps separate from rooms where gel electrophoresis of amplicons is performed.
- b) Use filter tips for micropipettes.
- c) Wear disposable powder-free gloves.
- d) Never store DNAnimal Kits and materials for DNA extraction together with samples or amplicons.
- e) Always perform extraction controls and PCR controls (NTCs).

12 / 32

🛟 eurofins

GeneScan

The following reactions are required for a RT-PCR run with 1, 8, 23 or 46 samples (incl. extraction controls). Add approx. 5% of volume as pipetting error compensation or use complete tubes.

Samples	1	8	23	46
NTCs	2	2	2	2
Pos. controls	2	2	2	2
Samples (duplicate)	2	16	46	92
Total no. reactions	6	20	50	96
Total MM volume	120 µL	400 µL	1 mL	2 mL
BasicMix	75 µL	250 µL	625 µL	1250 µL
			or	or
			1 tube	2 tubes
OligoMix	45 µL	150 µL	375 µL	750 µL
-			or	or
			1 tube	2 tubes
ROX (only for HR!)	2.7 µL	9 µL	22.5 µL	43.2 μL

7.2 Preparation of MasterMix

Calculate the number of reactions and amount of equivalent MasterMix volume before thawing and mixing the reagents and starting the practical work. The MasterMix is made of pre-made reagents ready for direct use after mixing of the two components.

Please be aware that the NTCs and positive controls (C+) are indispensable for the evaluation of the test and must be included in every run. PCR is performed in a volume of 25 μ L in 0.2 mL reaction tubes/plates according to the RT-PCR cycler instructions.

20 µL MasterMix for low ROX (LR) cyclers consist of 12.5 µL BasicMix + 7.5 µL OligoMix per reaction.

If you use one of the following cyclers, add also more ROXTMdye: ABI 5700, 7000, 7300, 7700, 7900, Opticon 2 or other cyclers laid out for high ROX (HR) concentration (please enquire in case of doubt). ROXTM is added increasing the volume of the MasterMix to 20.45 μ L.

20.45 μL HR MasterMix consist of 12.5 μL BasicMix + 0.45 μL ROX (50x) + 7.5 μL OligoMix per reaction.

13 / 32

DNAnimal Ident Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

- 1. Thaw only the volumes of BasicMix and OligoMix needed for analysis. Shake thoroughly.
- 2. Remove the required volumes and mix in a fresh tube. Freeze the rest.
- If the cycler requires a high ROX[™] level, add ROX[™].
- 4. Mix thoroughly.
- 5. The composed MasterMix can be stored up to 4 h in the refrigerator, but must not be used longer.

7.3 PCR Setup

Every run requires two no template controls (NTC) and two positive controls (C+). It is highly recommended to test the samples in duplicate.

- 1. Label all PCR reaction tubes.
- 2. Add 20 μL of the composed MasterMix to the wells.
- 3. Add 5 μ L of stabilisation buffer to NTCs.
- 4. Add 5 μL of control (sheep) DNA to positive controls.
- 5. Add 5 μL of sample DNA to test reactions.

NTC NTC

11a |11b

17a |17b

23a 23b

29a 29b

35a 35b

H 41a 41b 42a 42b

Plate layout for 46 samples.

NTC = no template control; C+ = positive control;

a and b = sample duplicates

B 5a 5b

С

D

G

C+

6a 6b

12a 12b

18a 18b

24a 24b

30a 30b

36a 36b

C+

1a 1b 2a

7a 7b

13a

19a |19b

25a

31a 31b

37a

2b 3a

8b 9a

14a 14b 15a 15b 16a

20a 20b 21a 21b 22a

26a 26b 27a 27b 28a

32a 32b 33a 33b 34a

43a | 43b | 44a | 44b | 45a | 45b | 46a | 46b

38a 38b 39a 39b 40a 40b

8a

13b

25b

37b

3b 4a 4b

9b 10a

GeneScan

DNAnimal Ident Sheep IPC (LR/HR)

Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

12

10b

16b

22b

28b

34b

7.4 Programming of Plate Documents

Before starting the practical work, program the plate document und the cycling conditions.

For description of the instrument programming please refer to the user manual of the respective instrument and software version.

Plate template

For the plate template, you may use our templates (send an e-mail mentioning your cycler model to kits@eurofins.com to receive a copy) or program your own template with the settings as below.

Cycling Conditions

The run parameters are to be programmed as follows:

Thermal Cycler Times and Temperatures			
1 HOLD	CYCLE (45 repeats)		
enzyme act.	denaturation	ann. & extension	
10 min at 95°C	15 sec at 95°C	90 sec at 60°C	
no data collection	no data coll.	data collection	

For other thermocyclers than the mentioned ones, it may be necessary to optimise the PCR parameters.

17 / 32

16 / 32

🛟 eurofins

GeneScan

DNAnimal Ident Sheep IPC (LR/HR)

Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

ABI 7500 (7500 Fast) / ABI 7900 HT

System	Detector	Reporter	Quencher
Sheep	8040a	FAM™	NONE
IPC	8040i	VIC™	NONE
Passive R	eference: RO	Х	

ABI 7500 Fast PCR is performed in the "Standard 7500" run mode.

ABI 7900HT Activate 9600 emulation mode

Bio-Rad CFX96 Touch™

System	Target Name	Dyes
Sheep	8040a	FAM™
IPC	8040i	VIC™

Passive Reference: not required

Ramp rate: default maximum ramp rate A reference dye is not used.

For further setup and evaluation information for your individual cycler, please see our application notes for your cycler model on our website www.eurofins.com/kits.

If you intend to use the Eurofins GeneScan evaluation sheet, please use the detector/assay names given below:

Stratagene Mx3005P (3000P)

System	Assay	Filter
Sheep	8040a	FAM™/SYBR® Green I
IPC	8040i	HEX™
Reference	ROX	ROX™

The fluorescence signals are scanned in the following order of filter: (1) ROX, (2) FAM, (3) HEX. The filter set gain settings for ROX, FAM and HEX are: pmt = 1x

<u>AriaMx</u>

System	Target Name	Dye
Sheep	8040a	FAM™
IPC	8040i	HEX™
Passive R	eference: ROX	

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

8. RESULTS

8.1 Evaluation

Refer to your cycler's manual for details. An evaluation (Excel[™]) sheet can be requested at kits@eurofins.com.

Threshold:

<u>ABI 7900HT, ABI 7500 (FAST), Mx3000P/Mx3500P:</u> The threshold should be placed in the region of exponential amplification across all of the amplification plots. This region is depicted in the log view of the amplification plots as the linear portion of the plot. The threshold line should neither be placed in the plateau phase nor in the initial linear phase of amplification. Thresholds are typically in the range 0.03 - 0.1.

<u>Bio-Rad CFX96 Touch:</u> Auto calculated threshold can be used when appropriate.

<u>Agilent AriaMx:</u> Auto calculated threshold with default Background based threshold settings can be used when appropriate. Cycle range: 5 thru 9

Sigma multiplier: 10

21 / 32

Roche LightCycler®480 Instrument I

Detection Format: Multi Color Hydrolyses Probe

System	Filter
Sheep	FAM (483-533)
IPC	VIC/HEX/Yellow555 (523-568)

Passive Reference: not required

Color Compensation for FAM and HEX is performed according to the manual.

FAM signal is collected using the FAM detector R6G signal is collected using the VIC/HEX /Yellow555 detector.

Ramp rate is set to 4.4 for heating and 2.2 for cooling. A reference dye is not used.

On LC480, Color Compensation has to be performed before the first test is run. Please use the Eurofins Color Compensation kit with cat. no. 5427200201.

Save the created file/document.

20 / 32

🛟 eurofins

GeneScan

DNAnimal Ident

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

8.2 Interpretation of Results

Export Ct values to the Excel[™] sheet provided (please request to kits@eurofins.com), or create an own evaluation sheet following the parameters below, where you will find an overview of the algorithms applied for scoring of the results in our Excel[™] Evaluation Sheet.

Definitions

Ct Cut-offs	sheep: Mean Ct (C+) +8
	IPC: Mean Ct (NTC) +3
Please Note:	For LightCycler 480 all amplification
For LC480:	curves with CP > 40 have to be checked
CP Cut-offs	visually for amplification
dR(n) Limits	sheep: Mean dR(n) (C+) x 0.2
	IPC: Mean dR(n) (NTC) x 0.33
(Mx3005P,	
Mx3000P,	
AriaMx,	For CFX Touch: End RFU
ABI 7500,	For LC480: Endpoint fluorescence
ABI7900)	
Outliers	Maximum acceptable outliers (C+): 0
	Maximum acceptable outliers (NTC): 0

Roche LC 480: No threshold settings

Baseline:

ABI RT-cyclers:	manual, 3-15 or automatic baseline
Stratagene Mx:	Adaptive
AriaMx:	Auto baseline correction can be
	used
Bio-Rad CFX96:	Baseline Subtracted Curve Fit
	Apply fluorescence drift correction
LC480:	no baseline setting necessary

Analysis mode: Roche LC 480

Apply the color compensation object (created for this analysis) to the experiment. To obtain CP values: use the "Abs Quant/2nd Derivative Max" analysis mode and High Sensitivity settings. To obtain endpoint fluorescence values: use the "Endpoint Genotyping" analysis mode.

DNAnimal Ident

Test reaction (sheep)

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

Inhibition control, scoring of IPC

Ct (CP for LC480) and dR(n)	Results
Ct_{IPC} sample $\leq Ct_{IPC}$ Cut-off and dR(n)_{IPC} sample $\geq dR(n)_{IPC}$ Limit	Sample valid
Ct_{IPC} sample > Ct_{IPC} Cut-off <u>or</u> dR(n) _{IPC} sample < dR(n) _{IPC} Limit	Sample invalid
No Ct _{IPC}	Sample invalid
Ct _{IPC} sample < Mean Ct _{IPC} NTC -3	Sample invalid

Ct	dR(n)	Result	
Ct sample ≤ Ct Cut-off	dR(n) sample ≥ dR(n) Limit	positive	
Ct sample ≤ Ct Cut-off	dR(n) sample < dR(n) Limit	Check amplification!	
Ct sample > Ct Cut-off	dR(n) sample ≥ dR(n) Limit	positive	
Ct sample > Ct Cut-off	dR(n) sample < dR(n) Limit	negative	
For LC480: Ct sample < Mean Ct (C+) -5	EF sample ≥ EF Limit	Check amplification!	
No Ct	-	negative	
Note: In case of "Check amplification!" the amplification			

plots must be checked visually for presence of a sigmoid PCR amplification signal.

24 / 32

🛟 eurofins

GeneScan

DNAnimal Ident

Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

8.3 Ambiguous results

If independently extracted DNAs show deviations in the results for "A" and "B", this may be due to sample material non-homogeneity or to non-uniformities in the DNA-extraction efficiency. Repeat DNA extraction and homogenize sample more thoroughly. If the ambiguous result persists, it is most likely due to a concentration close to the LOD.

Final result from combination of inhibition control and test reaction

IPC	Sheep test	Final result
Sample valid	Reaction positive	Positive
Sample valid	Reaction negative	Negative
Sample invalid	Reaction positive	Positive
Sample invalid	Reaction negative	Inhibited

Evaluation of the IPC

Calculate the MEAN Ct (Cp for LC480) Value from NTC. Refer to data from IPC-detector. To calculate the Ct cutoff, add 3 Ct.

Calculate the MEAN dR(n) Value from NTC. Refer to data from IPC-detector. The dR(n) cut-off f is 33% of the MEAN dR(n).

Evaluation of sheep specific test

Calculate the MEAN Ct Value from positive control. To calculate the Ct cut-off, add 8 Ct. Calculate the MEAN dR(n) value from positive control. The dR(n) cut-off is 20% of the MEAN dR(n).

Combine results of IPC and the sheep test to the final result.

25 / 32

DNAnimal Ident Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

8.4 Limit of detection

The relative limit of detection has been validated at 0.01% of sheep DNA in other species (total DNA content of 100 ng/rxn).

The absolute detection limit of the method is \leq 10 copies per reaction.

The limit of detection for the analysis of individual samples depends on several factors: the presence of PCR-inhibiting substances, the capability of the DNA extraction method to remove these inhibitors and the degree of DNA damage, and thus strongly on the type of sample and the DNA extraction procedure. Quantification is not possible with this kit.

9. PRODUCT USE LIMITATIONS

The GeneScan DNAnimal Kit is developed, designed, and sold for research purposes only. It is not to be used for diagnostic purposes or analysis of food and feed unless expressly cleared for that purpose by the appropriate regulatory authorities in the country of use. All due care and attention should be exercised in the handling of many of the materials described in this text.

28 / 32

🛟 eurofins

GeneScan

10. PRODUCT WARRANTIES AND SATISFACTION GUARANTEE

Eurofins GeneScan Technologies GmbH ("GeneScan") warrants the products manufactured by it will be free of defects in materials and workmanship when used in accordance with the applicable instructions and before the expiration date marked on the product packaging under the storage conditions, recommended in the instructions and/or on the package. Application protocols published by GeneScan are intended to be only guidelines for the buyers of the products. Each buyer is expected to validate the applicability of each application protocol to his individual application. GeneScan makes no other warranty, expressed or implied. There is no warranty of merchantability or fitness for a particular purpose.

GeneScan's sole obligation with the respect to the foregoing warranties shall be, at its option, to either replace or to refund the purchase price of the product(s) or part thereof that proves defective in materials or workmanship within the warranty period, provided the customer notifies GeneScan promptly of any such defect. GeneScan shall not be liable for any direct, indirect or consequential damages resulting from

29 / 32

DNAnimal Ident Sheep IPC (LR/HR) Test kit for qualitative detection of sheep DNA

Cat. No. 5422211710

12. TROUBLESHOOTING

Result	Possible mistakes/reasons	Possible verification and measures
No PCR signals from samples	Inhibition of PCR by inhibitory substances.	Clean DNA further* or dilute DNA solution.
	Inhibition by too much DNA.	Check DNA concentration/dilution.
No PCR signals from positive controls	Wrong PCR program.	Check and correct PCR program.
No amplification, neither from control DNA nor	MasterMix not properly prepared	Prepare fresh MasterMix, repeat PCR.
from sample DNA Positive PCR	Wrong PCR program. Contamination with	Check program. Optimise your
result for NTC	DNA/amplicons when mixing the PCR components.	precautions. Check your solutions. Decontaminate your equipment. Repeat the PCR.
Positive PCR result for extraction control	Contamination with sample material/DNA/	Check your solutions. Repeat extraction and
	amplicons/ during DNA extraction or PCR setup.	PCR.

* Repeat extraction of DNA from the sample. Repeat washing with 75 % ethanol three times. If necessary, clean DNA further e.g. with our DNA Cleaning Columns (cat. no. 5224700310).

economic loss or property damages sustained by buyer or any customer from the use of the product(s). A copy of Eurofins GeneScan Technologies GmbH terms and conditions can be obtained on request and is provided in our price lists.

11. IMPORTANT NOTES

- The TaqMan processes are covered by patents issued to Hoffmann-La Roche AG and Applied Biosystems Inc., which are applicable in certain countries. Eurofins GeneScan Technologies does not encourage or support the unauthorized or unlicensed use of these processes. Use of this kit is recommended for persons that either have a license or are not required to obtain a license.
- Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.